

Immersion Oil Compatibility With the ibidi Labware

As of July 25, 2024



ibidi Glass Coverslip Bottom Labware

When using ibidi Glass Bottom products with oil immersion objectives, there is no known incompatibility with any immersion oil on the market. All types of immersion oils can be used.

ibidi Polymer Coverslip (ibiTreat, Uncoated, and Protein-Coated) Bottom Labware

When using oil immersion objectives with the ibidi Polymer Coverslip, use only the immersion oils specified in the table below. The use of any non-recommended oil could damage the ibidi Polymer Coverslip, and the resulting leakage may harm objectives and microscope components. All immersion oils that are not listed in the table below should be considered non-compatible.

Company	Product	Ordering Number	Lot Number	Test Date
ibidi	ibidi Immersion Oil 2	50102	24-07-04	07/2024
Cargille	Type A	16482	100592	01/2017
Cargille	Type HF	16245	92192	01/2017
Carl Roth	Immersion oil	X899.1	414220338	01/2017
Leica	Immersion Liquid	11513859	n.a.	03/2023
Leica	Immersion Liquid Type G	11513910	n.a.	04/2024
Nikon	Immersion Oil F2 30cc*	MXA22192	n.a.	01/2020
Nikon	Silicone Immersion Oil 30cc	MXA22179	20191101	01/2020
Olympus	Silicone Immersion Oil	SIL300CS-30CC	N4190800	01/2017
Zeiss	Immersionol 518 F	444960-0000	220211	03/2023
Zeiss	Immersionol 518 F (30°C)	444970-9010	220816	03/2023
Zeiss	Immersionol 518 F (37°C)	444970-9000	220302	03/2023
Zeiss	Immersionol W 2010	444969-0000	101122	04/2012
Zeiss	Immersion Sil 406	444971-9000	80730	03/2023
Zeiss	Immersionol G	462959-9901	211117	03/2023

* Not available in the United States of America

